Application technologies

Laser Scanning Confocal Microscopy (**LSM**), Fluorescence Resonance Energy Transfer (**FRET**), Fluorescence Recovery After Photobleaching (**FRAP**), Low intensity fluorescence, Immunohistology, High Content Analysis/Screening (**HCA/HCS**), Phototoxicity studies, Hypoxia/Hyperoxia studies, Metabolic highly active cells

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Konfokale Laser Scanning Mikroskopie (**LSM**), Fluorescence Resonance Energy Transfer (**FRET**), Fluorescence Recovery After Photobleaching (**FRAP**), Schwach fluoreszierende Farbstoffe/Proteine, Immunhistologie, High Content Analysis/Screening (**HCA/HCS**), Fototoxizität, Exakte Kontrolle über Hypoxie/Hyperoxie, Stoffwechselaktive Zellen

Imaging Plate FC
24 well / 96 well

Features & benefits

Imaging Plates FC are black multiwell plates compliant to the SBS (Society for Biomolecular Screening) standard with a bottom made from a 25 µm fluorocarbon (FC) film. The thin film is stretched around each single well to ensure high level planicity. The FC film is gaspermeable and enables the fast equilibration of gas partial pressure in the cellular microenvironment with the surrounding air. In addition the film shows significant light transmission already at a wavelength of 210 nm.

Imaging Plates FC can be used for high resolution live cell microscopy. The plates are also suitable for a wide range of fixation and staining protocols (formaldehyde, glutaraldehyde, alcohols, acetone). Acetone should not be applied for periods longer than 10 minutes since the plate body is made from polystyrene and the plate body can detach from the bottom. Imaging Plates CG can be used at temperatures between -20°C and 50°C. Cell adhesion, spreading and distribution are improved due to our well established plasma surface modification technologies.

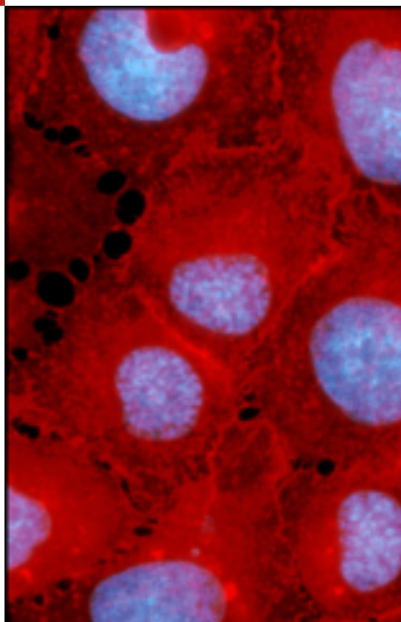
The thin FC film plate bottoms make the plates suitable for the application of high resolution immersion objectives. For objectives with magnifications $\geq 40\times$ objectives with a correction ring or collar have to be applied to get optimal focus control. Water, glycerine and oil can be used as immersion medium.

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Product highlights


- plate bottom **permeable for gases, UV-A and UV-B light**
- **high resolution** imaging for **throughput applications**
- 96 well plate with "**low-skirt design**", full bottom area can be used with immersion objectives
- **Advanced Tissue Culture** surface quality for cell adhesion and spreading

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

Fluorcarbon film properties:

Thickness / Stärke	25 µm
Thickness tolerance / Toleranz	± 10%
Flexible film is stretched around each single well	
Surface treatment / Oberfläche	Tissue Culture (TC)
Refractive index / Brechungsindex	1.34
Abbe's number / Abbe'sche Zahl	70
Light transmission @ 240nm	> 85%
Light transmission ≥ 300nm	> 90%
Oxygen permeability	>6300 cm ³ /(m ² *d* bar)
CO ₂ permeability	>7000 cm ³ /(m ² *d* bar)

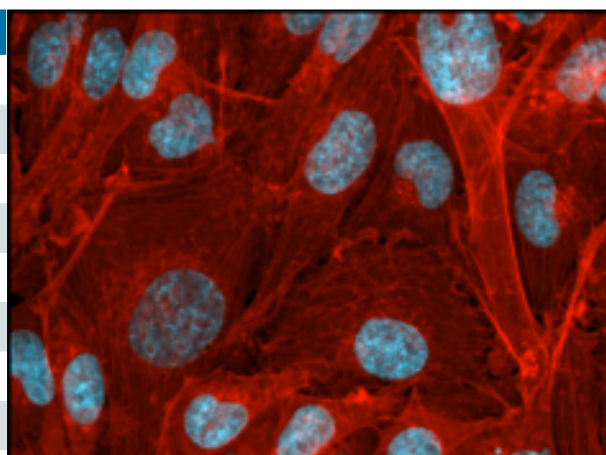Imaging Plate FC 24 well / 96 well

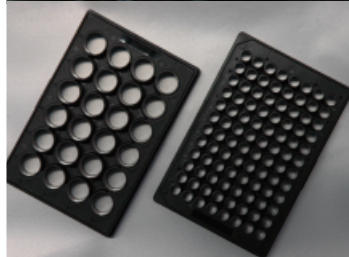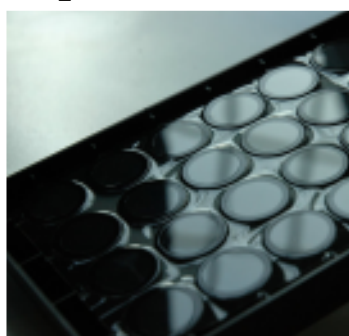Article number / Artikelnummer	3231	3241
Format	24 well	96 well
Plate length	127,76 mm	127,76 mm
Plate width	85,48 mm	85,48 mm
Inner well diameter (imaging plane)	13.2 mm	6 mm
Distance well to well center	18 mm	9 mm
Distance center well A1 plate border (long side)	15.74 mm	11.24 mm
Distance center well A1 plate border (small side)	18.88 mm	14.38 mm
Distance plate bottom to image focus plane	2.5 mm	400 µm
Inner well flatness	≤ 10 µm	≤ 10 µm
Well to well flatness	≤ 50 µm	≤ 50 µm
Total well volume	1880 µl	428 µl
Suggested working volume	500-1000 µl	100-200 µl

The polymer bodies of all labware react to temperature shifts with changes of their dimension. Please ensure that you can achieve equilibrated environmental temperature conditions during microscopy to avoid drift of the focal plane.

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General information

All Imaging Plates are shipped single packed with lid.
All lids provide condensation rings and spacers for optimal ventilation.

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zell-kontakt GmbH