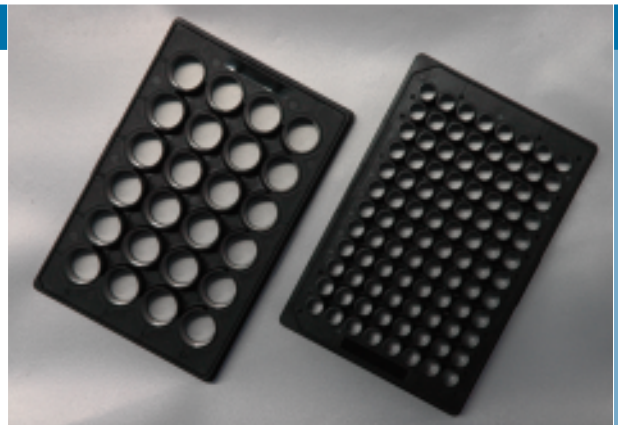
Application technologies

Laser Scanning Confocal Microscopy (LSM), Total Internal Reflection Fluorescence (TIRF), Differential Interference Contrast (DIC), Fluorescence Correlation Spectroscopy (FCS), Fluorescence Resonance Energy Transfer (FRET), Fluorescence Recovery After Photobleaching (FRAP), Low intensity fluorescence, Fluorescence In-Situ Hybridization (FISH), Immunohistology, High Content Analysis/Screening (HCA/HCS)

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Konfokale Laser Scanning Mikroskopie (LSM), Total Internal Reflection Fluorescence (TIRF), Differentieller Interferenz Kontrast (DIC/Nomarski), Fluoreszenz Korrelationsspektroskopie (FCS), Fluoreszenz Resonance Energy Transfer (FRET), Fluoreszenz Recovery After Photobleaching (FRAP), Schwach fluoreszierende Farbstoffe/Proteine, Fluoreszenz In-Situ Hybridisierung (FISH), Immunhistologie, High Content Analysis/Screening (HCA/HCS)

Imaging Plate CG
24 well / 96 well

Features & benefits

Imaging Plates CG are black multiwell plates compliant to the SBS (Society for Biomolecular Screening) standard with a bottom made from borosilicate cover glass. The plate bodies are made from polystyrene and are tightly sealed against the glass bottoms by adhesive.

Imaging Plates CG can be used for high resolution live cell microscopy. The plates are also suitable for a wide range of fixation and staining protocols (formaldehyde, glutaraldehyde, alcohols, acetone). Acetone should not be applied for periods longer than 10 minutes since the plate body is made from polystyrene and the plate body can detach from the glass bottom.

Imaging Plates CG can be used at temperatures between -20°C and 50°C.

Cell adhesion, spreading and distribution are improved due to our well established plasma surface modification technologies.

The cover glass plate bottoms make the plates suitable for the application of high resolution immersion objectives. Water, glycerine and oil can be used as immersion medium.

Imaging Plates CG are available in 24 well and 96 well format.

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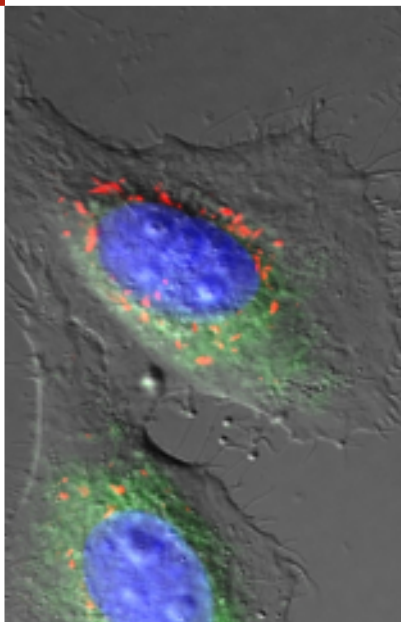
Product highlights


- **extraordinary planicity** of the imaging plane
- **high resolution** imaging for **throughput applications**
- 96 well plate with "**low-skirt design**", full bottom area can be used with immersion objectives
- **Advanced Tissue Culture** surface quality for cell adhesion and spreading

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

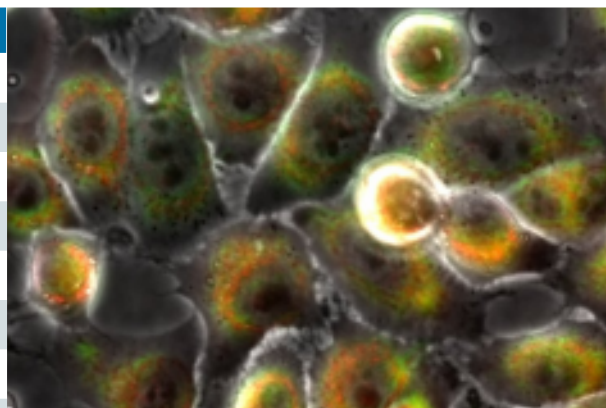
Imaging Plate CG 24 well / 96 well

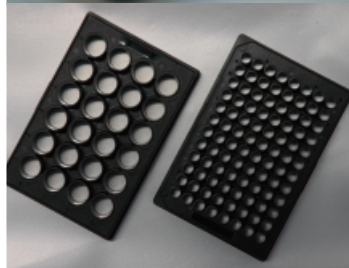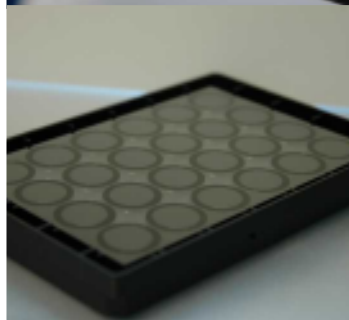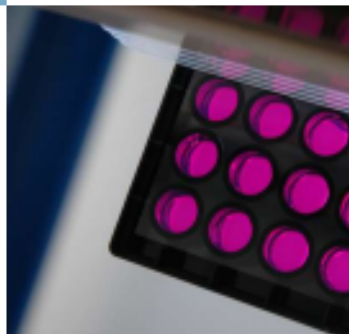Article number / Artikelnummer	5231	5241
Format	24 well	96 well
Plate length	127,76 mm	127,76 mm
Plate width	85,48 mm	85,48 mm
Inner well diameter (imaging plane)	13.2 mm	6 mm
Distance well to well center	18 mm	9 mm
Distance center well A1 plate border (long side)	15.74 mm	11.24 mm
Distance center well A1 plate border (small side)	18.88 mm	14.38 mm
Distance plate bottom to image focus plane	2.5 mm	400 µm
Inner well flatness	≤ 10 µm	≤ 10 µm
Well to well flatness	≤ 50 µm	≤ 50 µm
Total well volume	1880 µl	428 µl
Suggested working volume	500-1000 µl	100-200 µl

The polymer bodies of all labware react to temperature shifts with changes of their dimension. Please ensure that you can achieve equilibrated environmental temperature conditions during microscopy to avoid drift of the focal plane.

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General information

All Imaging Plates are shipped single packed with lid.
All lids provide condensation rings and spacers for optimal ventilation.

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**zell-kontakt GmbH**